

# I N S T R U C T I O N S

## Mercury Collector Replacement Pads

Catalog No. F36251-0000

These open cell polyurethane pads are replacements for the pad supplied with the Bel-Art mercury collector F32650-0000

When handling the used pad, wear gloves and a mask to prevent contact with hazardous mercury. Dispose of the used pad in accordance with local regulations.

1. Using a scraper, such as a putty knife, remove the used pad from the three legged platform and scrape off excess dried cement.
2. To install a new pad on the platform, use rubber cement (usually available at art supply stores). Apply a layer of cement to the platform and center the pad on the platform with light pressure. For your safety, comply with the warnings on the rubber cement instructions.

